



Advantech Manufacturing, Inc

2450 S. Commerce Drive
New Berlin, WI 53151

Toll Free: 800.511.2096

Tel: 262.786.1600

Fax: 262.786.5074

www.advantechmfg.com
sales@advantechmfg.com

CARE AND CLEANING FOR METAL FRAMED WIRE CLOTH SIEVES

*(For care and handling of electroformed nickel precision sieves,
please refer to: CARE AND CLEANING OF PRECISION SIEVES.)*

- Clean sieves from **BOTTOM** side with a brush. Use extreme care to avoid damage.
- Never use sharp objects to clean your wire cloth sieves.
- Wash in a mild solution of detergent and warm water. For a more thorough cleaning, some ultrasonic* cleaners may be used. (**Caution: Refer to your ultrasonic manufacturer instructions prior to use as some devices may cause screen damage.*)
- **CAUTION:** Do not expose sieves to temperatures above 261° Fahrenheit as solder will begin to soften.
- Do not use forced air to dry #100 or finer wire cloth mesh sieves. This may result in screen damage.
- Store equipment in a clean, dry environment when not in use.